



ECONOMICAL DISINFECTION SUITABLE FOR USE IN LOWER GRADE AREAS



Klercide™ 70|30 Filtered Alcohol

Cleanroom filtered 70|30 v/v alcohol, either IPA or Denatured Ethanol, blended with deionised water

Klercide Filtered Alcohol is available as either 70% v/v Isopropyl Alcohol blended with deionised water or 70% v/v Denatured Ethanol blended with deionised water.

The alcohol blend is 0.2 micron filtered, then filled and double bagged in a Grade C (ISO Class 7) cleanroom.

All spray bottles incorporate the patented SteriShield Delivery System and an adjustable trigger spray. The unique way in which the SteriShield Delivery System works, as a closed system, ensures that the integrity of the contents is preserved throughout use.

Available in 1 litre trigger spray and 5 litre capped bottles.

Klercide 70|30 Alcohol contains no propellants, so does not require any special disposal procedures.

The range:

- ▲ Includes trigger sprays, fitted with the SDS trigger spray, to give an in-use shelf life of 3 months
- ▲ Is available in a variety of formats
- ▲ Incorporates a fully adjustable trigger spray on the 1 litre bottle



Klercide™ 70|30 Filtered Alcohol



Cleanroom filtered 70|30 Filtered Alcohol v/v alcohol, either IPA or Denatured Ethanol, blended with deionised water

Why use Klercide 70|30 Filtered Alcohol?

EASY TO USE

Convenient range of easy to use products are available in 1 litre spray or 5 litre capped formats. The large droplet size of the spray reduces the risk of inhalation.

COST EFFECTIVE

All of the alcohol can be dispensed from the bottle so there is no wastage. The SteriShield Delivery System has been validated to give an in-use shelf life of three months, so there is no need to discard unused product.

OPTIMUM SURFACE COVERAGE

The fully adjustable trigger spray allows the liquid to be dispensed as either a jet or spray enabling thorough wetting of the surface so that disinfection procedures are effective.

PROTOCOL

Ready-to-use so can be sprayed directly onto the surface to be sanitised. Always ensure the whole surface is in contact with the liquid. For optimum results wipe the surface to physically remove any contaminants.

VALIDATION

Manufactured by an ISO 9001:2015 and ISO 14001:2015 certified company.

EN test data available on request.

Certificates of Analysis and Irradiation are available for every batch.

Product has a two year shelf life.

Full supporting documentation is available on this product.

**Effective.
Economical.
Efficient.**

APPLICATION AREA

For use on floors, walls and hard surfaces.

INSTRUCTIONS FOR USE

Spray: For optimum results hold bottle upright and spray from a distance of 10 cm to 20 cm. Always close the nozzle after use.

Wetting: For optimum results ensure complete wetting of the surfaces using the spray and wipe technique (max. 40 ml/m²).

INGREDIENTS

100 ml contains: Either 70% v/v Isopropyl Alcohol blended with deionised water or 70% Denatured Ethanol blended with deionised water.

PACKAGING INFORMATION

Klercide 70|30 Alcohol (Filtered)

UNIT SIZE	UNIT OF SALE
1 L spray	6 x 1 L
5 L capped	4 x 5 L

FOR FULL PRODUCT ORDER CODES, PLEASE VISIT OUR WEBSITE.



70|30 Denatured Ethanol
1 L SDS Spray, 5 L Capped



70|30 IPA
1 L SDS Spray, 5 L Capped

USE BIOCIDES SAFELY, ALWAYS READ THE LABEL AND PRODUCT INFORMATION BEFORE USE

WORLDWIDE HEADQUARTERS

1 Ecolab Place
St. Paul, MN 55102
USA

www.ecolab.com/lifesciences

EUROPE HEADQUARTERS

Richtistr. 7
8304 Wallisellen
Switzerland

www.ecolablifesciences.com

